

# Characterization of Surfaces that Generate Structural Color by Total Internal Reflection

Our research involves the design of microstructures that generate interference colors. Our first report of this optical mechanism appeared in Goodling, A. E., et al. (2019). "Colouration by total internal reflection and interference at microscale concave interfaces." *Nature* 566(7745): 523-527. Light bounces by total internal reflection within the microstructures and optical interference occurs between light rays undergoing different paths of reflection.

We design microstructured surfaces to control the reflection of the light and the colors produced. We need to measure the surface topography to be able to correlate optical properties with the surface geometries. The geometries of the microstructures, such as their depth, width, angles, and surface roughness impact the optical properties. We use the optical profilometer to characterize the surface profiles of the microstructures and the uniformity across the surface. Sometimes our structures have steep angles, so being able to use different modes of surface profiling is important to get the most accurate determination of the microstructure geometry.



[Penn State University,](#)  
[Department of Chemistry](#)




and Department of Materials Science and Engineering. The Department of Chemistry at Penn State offers excellence in the chemical sciences through a vibrant and collaborative academic community.

Authors: Lauren Zarzar, Nate Sturniolo, Caleb Meredith, and Amy Goodling.



**PennState**

■ Measurements

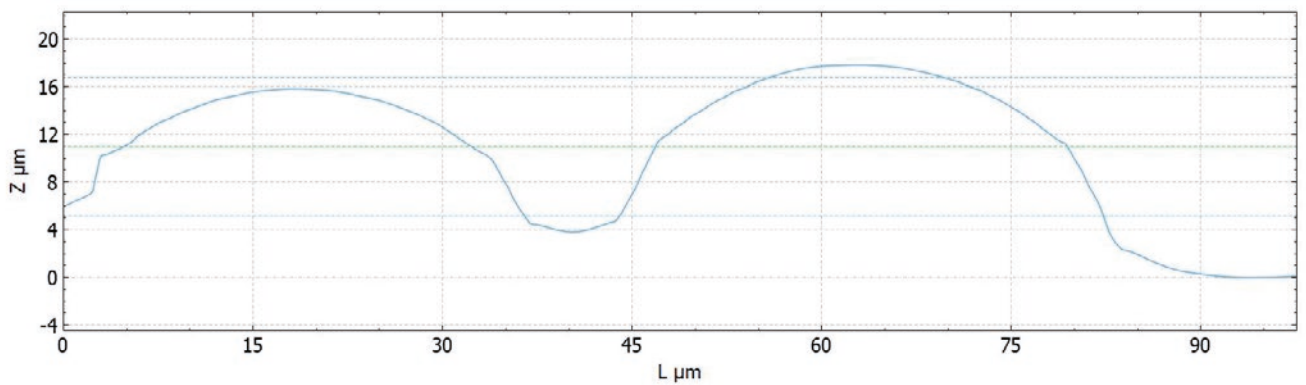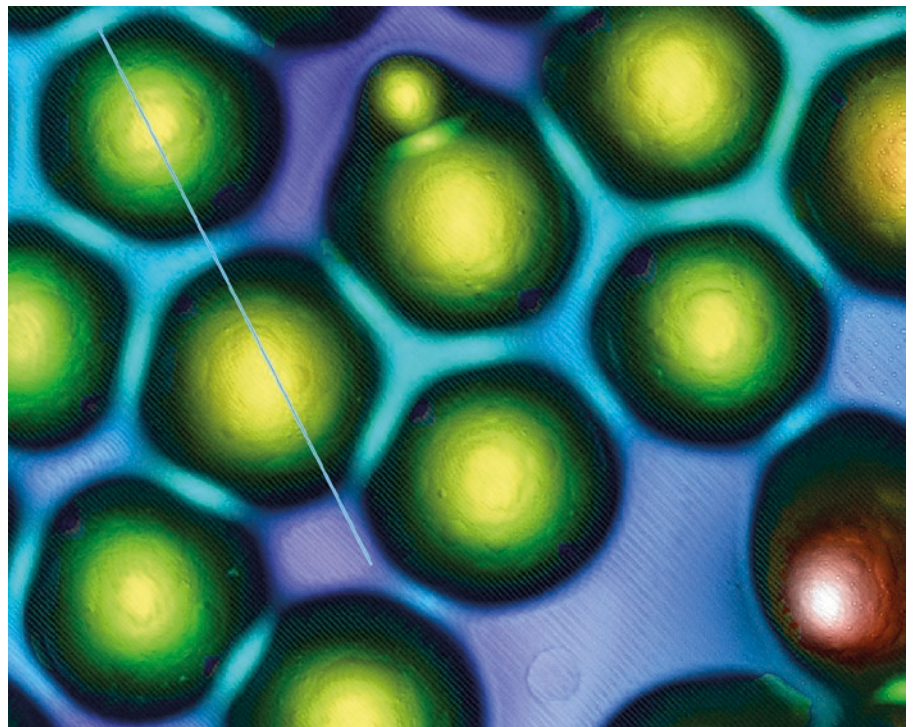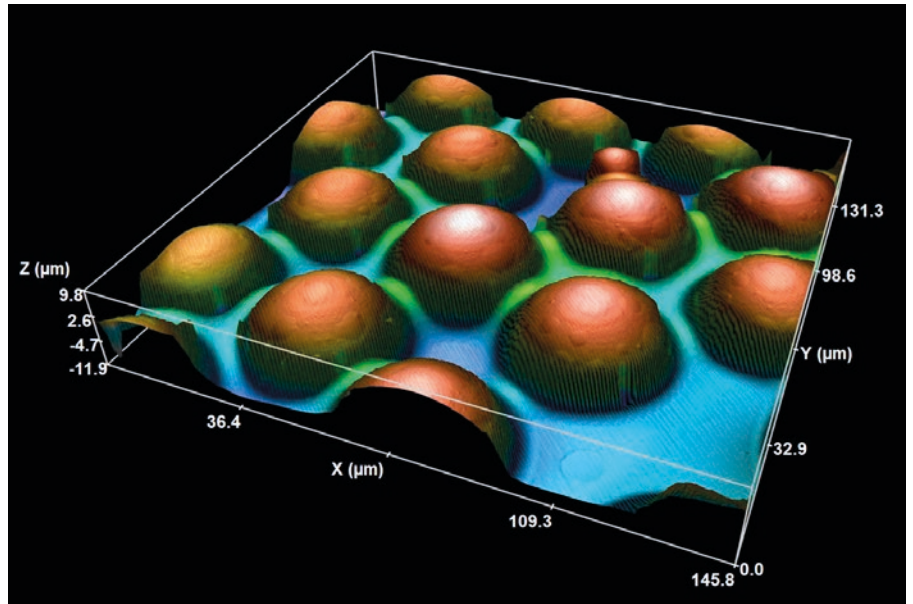

---

## ■ Conclusions

Using the profilometer, we have been able to image our microstructured surfaces with high spatial resolution in XYZ to yield important information regarding the surface roughness of our samples, reproducibility after structural replication, and variation across the surface that may impact the optical properties. Sensofar's lynx profilometer was the best option for us because of the versatility in imaging modes available. Sometimes we need high accuracy in the height distribution across the surface and choose to use white light interference. Other times we need to get information on the sidewalls of the microstructures that have relatively steep angles, so confocal imaging is better. By choosing to use Sensofar's lynx, we have access to both of these imaging modalities which means we can get even more information about our sample than just using a confocal profiler or white light interferometer on its own. Sensofar worked closely with us to choose the specific profilometer and add-on options that worked best for specific samples and imaging needs.

## ■ References

1. Goodling, A. E., et al. (2019). **Colouration by total internal reflection and interference at microscale concave interfaces**. *Nature* 566(7745): 523-527.
2. Goodling, A. E., et al. (2020). **Tunable and Responsive Structural Color from Polymeric Microstructured Surfaces Enabled by Interference of Totally Internally Reflected Light**. *ACS Materials Letters* 2(7): 754-763.



SENSOFAR is a leading-edge technology company that has the highest quality standards within the field of surface metrology

Sensofar Metrology provides high-accuracy optical profilers based on confocal, interferometry and focus variation techniques, from standard setups for R&D and quality inspection laboratories to complete non-contact metrology solutions for in-line production processes. The Sensofar Group has its headquarters in Barcelona, known as a technology and innovation hub in Europe. The Group is represented in over 30 countries through a global network of partners and has its own offices in Asia, Germany and the United States.

**HEADQUARTERS**

**SENSOFAR METROLOGY** | BARCELONA (Spain) | T. +34 93 700 14 92 | [info@sensofar.com](mailto:info@sensofar.com)

**SALES OFFICES**

**SENSOFAR ASIA** | SHANGHAI (China) | T. +86 021 51602735 | [info.asia@sensofar.com](mailto:info.asia@sensofar.com)

**SENSOFAR GERMANY** | MUNICH (Germany) | T. +49 151 14304168 | [info.germany@sensofar.com](mailto:info.germany@sensofar.com)

**SENSOFAR USA** | NEWINGTON (USA) | T. +1 617 678 4185 | [info.usa@sensofar.com](mailto:info.usa@sensofar.com)

[sensofar.com](http://sensofar.com)